DNA from fruit


Komal Zahrah, Henriette Mair-Meijers, Justyna Resztak, Anthony Findley, Francesca Luca

Center for Molecular Medicine and Genetics
Wayne State University
fluca@wayne.edu

DNA from Fruit

Materials

- 1 Resealable (e.g. Ziplock) bag
- 1 Banana or 2 strawberries (can be substituted by a tomato, kiwi, apple)
- 2 tsp Dish detergent
- ½ Cup of water
- 2 Clear cups (plastic or glass)
- 1 Coffee filter or fine mesh strainer
- 1 tsp of salt
- ½ Cup of cold rubbing alcohol (Put bottle in freezer before use)
- Coffee stirrer or wooden skewer
- Measuring spoons

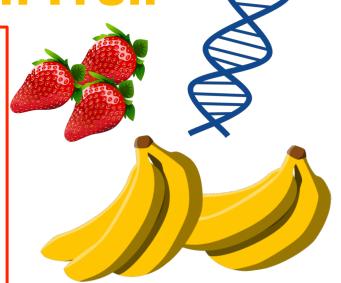

Procedure


- 1. Take the peel off the banana (or use 2 strawberries).
- 2. Put the fruit into the plastic bag.
- 3. Seal the bag.
- 4. Completely mash the fruit with your hands (this breaks open the cells and releases the DNA!).
- 5. In a cup, make the DNA extraction liquid:
 - 1. Mix together 2 teaspoons of detergent, 1 teaspoon of salt, and ½ cup of water
- 6. Add 2 teaspoons of the DNA extraction liquid into the bag with the fruit.
- 7. Reseal the bag and gently smash/mix for another minute (avoid making soap bubbles).
- 8. Place the coffee filter/strainer inside the other cup.
- 9. Open the bag and pour the content from the bag into the filter.
- 10. Gently squeeze the filter to let the liquid seep through.
- 11. Next, pour down the side of the cup an equal amount of cold rubbing alcohol as there is fruit liquid and do not mix or stir.
- 12. Within a few seconds, watch for the development of a white cloudy substance (DNA) in the top layer above the fruit extract layer.
- 13. Tilt the cup and pick up the DNA using a plastic coffee stirrer or wooden stick.
- 14. You have now isolated DNA from fruit!

What is DNA?

"It's a history book - a narrative of the journey of our species through time

It's a shop manual, with an incredibly detailed blueprint for building every human cell.

And it's a transformative textbook of medicine, with insights that will give health care providers immense new powers to treat, prevent and cure disease." - Francis Collins

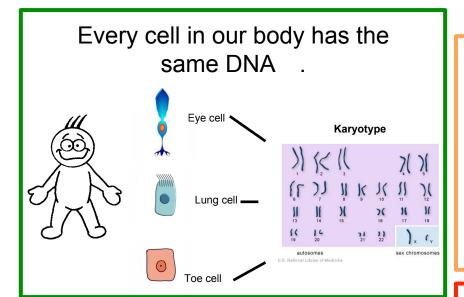

WHAT CAN WE LEARN FROM THE HUMAN GENOME?




Better understanding of human disease

Personalized medicine & Pharmacogenetics






Greater insight into cognitive function

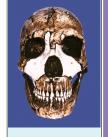

Insight into human origins



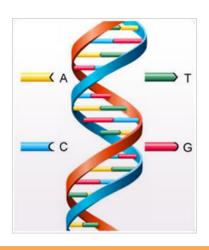


How much DNA is in one human cell?

Genome = 46 (23 x 2) chromosomes Genome = approx. 3 billion

base pairs. If stretched = 2 meters (6.6 ft) DNA sequence in any two people is 99.9% identical – only 0.1% is unique!

How much DNA is in one strawberry cell?

Genome = 56 (7 x 8) chromosomes Genome = approx. 960 million base pairs

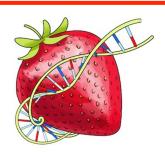

"Silly Science" is organized by the Luca Lab.



www.lucalab.wayne.edu School of Medicine



Credit

- Banana picture:

 https://upload.wikimedia.org/
 wikipedia/commons/thumb/f/f7/

 Bananas.svg/560px-Bananas.svg.png
- Strawberry picture:
- https://pngimg.com/uploads/ strawberry/strawberry PNG2586.png
- Dna picture: <u>https://pngimg.com/uploads/dna/dna PNG15.png</u>
- Adapted from:

 https://www.genome.gov/Pages/
 Education/Modules/

 StrawberryExtractionInstructions.pdf